



SOLO-X update procedure.

- Close all applications and disconnect the SOLO-X from the PC.
Insert the update mode connector.




- Install the latest USB drivers:
http://www.gentec-eo.com/download/software-and-firmware/usb-drivers/Monitor_USB_DRIVERS.exe
- Power-OFF the SOLO-X if it is ON.
- Connect the SOLO-X to the PC USB port using the cable.
- Open the device manager:

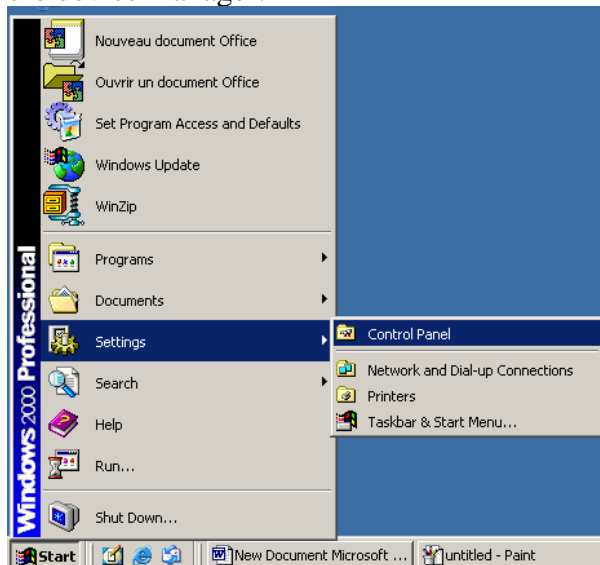

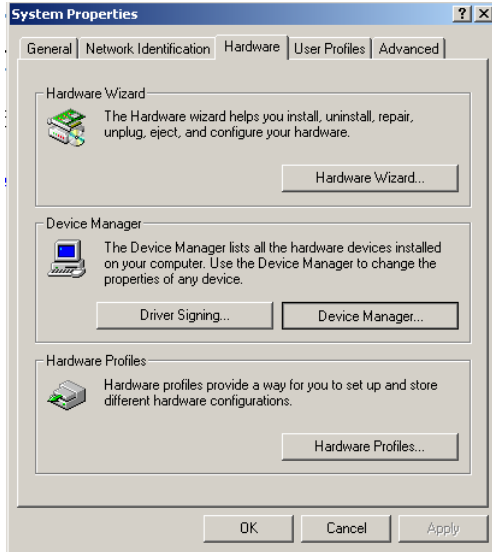

- Note which COM port is used by « Prolific USB-to-Serial» (com 3 in this case). The current Firmware Updater supports only COM ports under 9.

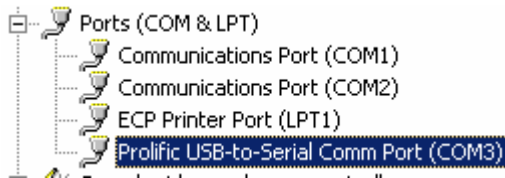

- Execute the SOLO update software:
<http://www.gentec-eo.com/download/software-and-firmware/firmware/Solo-V3.00.08.exe>

- Select the COM port found above and click on "Test communication with device". If successful, click on "Start Update":



- When update is over, close the Firmware Updater.
- Remove the update connector.
- Reboot the SOLO-X by turning it OFF and back ON.
- Install the latest version of PC-SOLO:

<http://www.gentec-eo.com/download/software-and-firmware/software/PC-Solo-V3.02.04.exe>